



UNPACKING

Carefully remove your Scout balance and each of its components from the package. Store the packaging material to ensure safe storage and future transport.

Included Components:

- Balance
- User Guide & Instructions
- Warranty Card
- AC Adapter & Plug
- Pan
- Calibration Mass (only for SJX/E models with 64g capacity)
- Scoop (only for models with 64g capacity)

CALIBRATION

When the balance is operated for the first time, a calibration is recommended to ensure accurate weighing results.

• External Span Calibration

With balance turned on:

Step-1	Press and hold Menu until [MENU] (Menu) is displayed.
Step-2	Release the button, the display will show [CAL].
Step-3	Press Yes to accept. [SPAN] will then be shown.
Step-4	Press Yes to begin the span calibration. Zero reading is stored automatically.
Step-5	Place the specified span calibration mass on the pan. Then, press Yes to accept the weight or No to select an alternate weight. [--C--] will be displayed while the reading is stored. The display will show [done] if the calibration was successful. The balance returns to the previous application mode and is ready for use.

• Internal Calibration (not available for SJX/E models)

With balance turned on:

Step-1	Press and hold Menu until [MENU] (Menu) is displayed.	
Step-2	Release the button, the display will show [CAL].	
Step-3	Press Yes to accept. [InCAL] will then be shown.	
Step-4	Press Yes to begin the Internal calibration. Zero reading is stored automatically.	
Step-5	Pull the InCal handle while the display show [PULL].	
Step-6	Push the InCal handle while the display show [PUSH].	
Step-7	Repeat the step-5, and step-6 one more time.	
Step-8	The display will show [done] if the calibration was successful. The balance returns to the previous application mode and is ready for use.	

OPERATION & APPLICATIONS

The balance can be configured to operate in one or more Application modes. Press and hold **Mode** key to select your Application mode.

The Scout SJX balance incorporates the following Applications:

[WEIGH]	[COUNT]	[PERCENT]	[CHECK]	[TOTAL]	[SPEC.Gr]	[Hold]
Weighing	Parts Counting	Percent Weighing	Check Weighing	Totalization	Density Determination	Display Hold

ADDITIONAL FEATURES


• **Connecting the Interface**

Use an optional interface connectivity kit to connect the balance either to a computer, printer or OHAUS auxiliary display.

Below Interface kit accessories are available:

RS232, USB Host, USB Device, Ethernet, Bluetooth.



Interface connection on the rear of the balance:

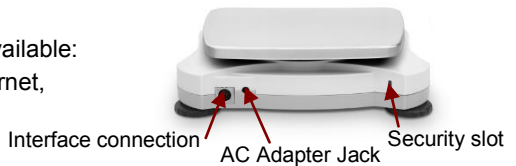

* Interface kits may vary according to local regulations

• **Weigh Below Hook**

The weigh below hook is located at the reverse side of the battery cover as shown. To use this feature, remove the red protective cover underneath for the weigh below opening. For SJX, the weigh below hook is integrated and located under the protective cover. For SJX/E, the weigh below hook is stored in the battery compartment (back side of the battery cover).

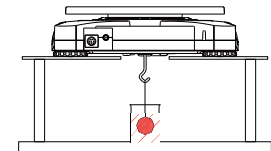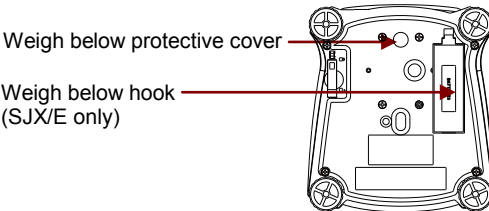

INSTRUCTION MANUAL

For more detailed information about the balance, please refer to the product instruction manual. The instruction manual can be obtained either by scanning the QR code or by going to the Scout SJX product page to download it.

Go to www.ohaus.com to register your product's warranty.

Product Page: <http://www.ohaus.com/>

QR Code:

