

Visit www.scaleman.com to see more Mettler-Toledo Balances we offer

Menu Map

Menu Operation

- Enter menu
- Leave menu with saving
- Enter menu selection
- Scroll down
- Scroll up
- One step back
- Leave menu without saving

Main Menu

- BASIC → Basic Menu
- ADVANCE → Advanced Menu
- INT.FACE → Interface Menu
- PROTECT → Protection

Protection

Menu	Selection
PROTECT	OFF
	ON

Basic Menu

Topic	Selection
DATE	01.01.11
TIME	+1H
	-1H
SET.TIME	12:00
1:10	OFF
	ON
UNIT 1	g
	kg
UNIT 2	g
	mg
SET ID	
PRT.MENU	
RESET	NO?
	YES?

Advanced Menu

Topic	Selection
ENVIRON	STD
	UNSTAB
	STABLE
CAL	ADJ.OFF
	ADJ.SET → 200.00 g
DATE.FRM	DD.MM.Y
	MM.DD.Y
	Y-MM-DD
	DDMM.Y
	MMDD.Y
TIME.FRM	24.MM
	12.MM
	24.MM
	12.MM
RECALL	OFF
	ON
STANDBY	RON → 60
	ROFF
B.LIGHT	B.L ON
	B.L OFF
AZERO	ON
	OFF
SRV:ICON	OFF
	ON
SRV:RST	NO?
	YES?

Interface Menu

Topic	Selection
R5232	PRINTER → PRT.STAB
	PRT.AUTO
	PRT.ALL
PC.DIR	PRT.STAB
	PRT.AUTO
	PRT.ALL
HOST	SNJ.OFF
	SNJ.STJ
	SNJ.CONT
	SNJ.AUTO
	SNJ.ALL
	2.DISP
HEADER	NO
	DATE.TM
	DATE.BAL
SINGLE	NET
	G/TM
SIGNAL	OFF
	ON
LN.FEED	DD
ZERO.PRT	OFF
	ON
COMSET	MT-SICS
	SAPT
BAUD	9600
	4800
	600...38400
BIT.PAR	8:NO
	7:NO
	7:MARK
	7:SPACE
	7:EVEN
	7:ODD
STOPBIT	1 BIT
	2 BITS
HD.SHK	ON:OFF
	P.TS.CTS
	OFF
P5.TE.O.L.	CR LF
	CR
	LF
	TAB
P5.CHAR	IBM.DOS
	ANSI.WIN
INTERVL	00000

Legend

- Factory setting
- Only those weight units allowed by the appropriate national legislation are selectable
- Not available with approved models
- Not available with approved models with e = d
- Settings are automatically set for the 2nd display
- Only visible if "PRINTER" is selected
- Only visible if "HOST" is selected
- Only visible if "PC-DIR" is selected
- Only visible if "PRT.AUTO" is selected

Visit www.balances.com to purchase your Mettler-Toledo Jewelry Scales